

TZM0880 Series 0.8-8.0X Monocular Zoom Objective (MZO)



Contents

| | |
|---|----|
| 1 The Basic Characteristic of TZM0880 Series MZO | 2 |
| 1.1 Introduction of TZM0880 Series MZO..... | 2 |
| 1.2 The Modular Design of TZM0880 Series MZO | 2 |
| 1.3 The Nomenclature of TZM0880 Series MZO | 3 |
| 1.3.1 Module One: Middle Zoom Module | 3 |
| 1.3.2 Module Two: Auxiliary Lens Module | 3 |
| 1.3.3 Module Three: TV Lens..... | 3 |
| 1.3.4 Module Four: Bracket Adapter | 4 |
| 1.3.5 Module Five: The Optional Module..... | 4 |
| 1.3.6 The Nomenclature of TZM0880 Series MZO | 6 |
| 1.3.7 The Configuration of TZM0880 with Five Modules..... | 7 |
| 1.4 The Basic Characteristics of TZM0880 Series MZO | 7 |
| 1.5 The Optical Specifications of TZM0880 with Different Auxiliary Lens and TV Lens | 8 |
| 1.6 The Dimension of TZM0880 Series MZO | 9 |
| 1.7 How to Configure TZM0880 Series MZO | 10 |
| 2 The Connection of TZM0880 Series MZO and Camera | 11 |
| 3 Pictures Taken with TZM0880 Series MZO | 11 |
| 4 TZM0880 Series MZO Packing List | 12 |
| 5 Order List of TZM0880 Series MZO | 14 |
| 6 Service | 15 |

1 The Basic Characteristic of TzM0880 Series MZO

1.1 Introduction of TzM0880 Series MZO

TzM0880 series MZO (Monocular Zoom Objective) is an ideal choice for machine vision, industrial inspection, scientific research and education. The design is based on the bilateral parallel light path principle and the modular design concept, with superior optical performance and strong compatibility.



Figure 1 TzM0880 with different module

1.2 The Modular Design of TzM0880 Series MZO

The modular design of **TzM0880** series MZO is showed in Figure 2. The design consists of five basic modules, including **Middle Zoom Module**, **Auxiliary Lens Module**, **TV Lens Module**, **Bracket Adapter Module** and **Optional Module**.

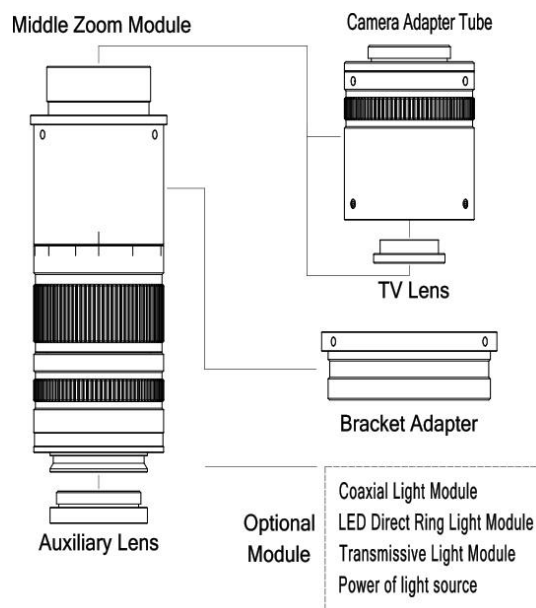


Figure 2 The five basic modules of TzM0880 series MZO.

1.3 The Nomenclature of TZM0880 Series MZO

Because of TZM0880 series MZO's modular design, each module of TZM0880 series MZO is named respectively. The specific naming rules and corresponding parameters of each module are as follows:

1.3.1 Module One: Middle Zoom Module

The Middle Zoom Module (code TZM0880a), which is the main body of TZM0880 series MZO. TZM0880 represents the code of Middle Zoom Module and letter a represents the auxiliary code of Middle Zoom Module. Details are listed in Table 1.

Table 1 The order number and function of the Middle Zoom Module

| Order Number | Meaning | TZM0880 | Postfix meaning |
|--------------|--|-----------------|-----------------|
| TZM0880 | Ordinary Middle Zoom Module | ZM: Zoom | NA |
| TZM0880D | Middle Zoom Module with detent to fix the magnification function | 0880: 0.8X~8.0X | D: Detent |

1.3.2 Module Two: Auxiliary Lens Module

The replaceable Auxiliary Lens (code W3bbb), which can be threaded to the bottom of the Middle Zoom Module, here 3 is the series number, bbb represents the magnification of Auxiliary Lens. Different magnification of Auxiliary Lens defines the entire magnification of TZM0880 series MZO, resulting different working distance as well. For example, W3100 Auxiliary Lens with 1.00x magnification reaches 86mm working distance, while W3050 Auxiliary Lens will have a reduced magnification to half and the working distance will be extended to about 174mm. Auxiliary Lens and its corresponding optical parameters for TZM0880 are shown in Table 2.

Table 2 The order number and optical parameters of Auxiliary Lens

| Order number | Magnification | Working distance (mmm) |
|--------------|---------------|------------------------|
| W3035 | 0.35X | 248 |
| W3050 | 0.50X | 174 |
| W3075 | 0.75X | 115 |
| W3100 | 1.00X | 86 |
| W3150 | 1.50X | 56 |

Note: The same Auxiliary Lens with different TV lens is just used to adapt to different size image sensor, which will not have much impact on the field of view of the MZO.

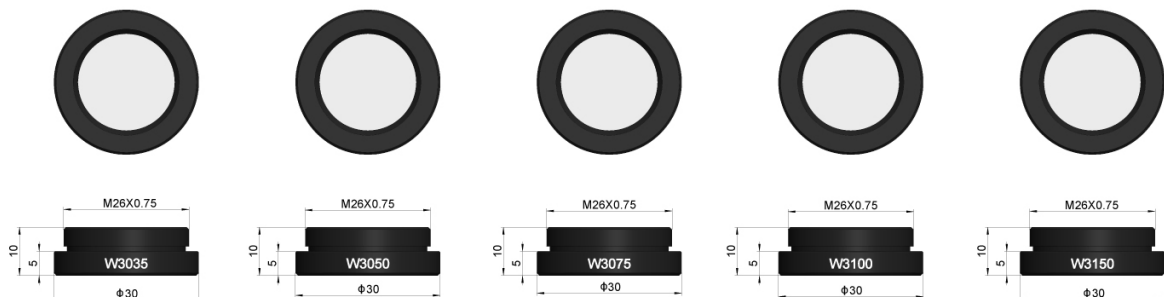


Figure 3 Auxiliary Lens for TZM0880 MZO

1.3.3 Module Three: TV Lens

The TV Lens (code TV3ccc), which can be threaded to the top of the Middle Zoom Module, here 3 is the series number, ccc is the magnification of the TV Lens. TV Lens with different magnification is used to adapt to sensors with different size. For example, the same field of view

TZM0880 Series MZO Help Manual
could be obtained when TV3100 is used with a 2/3" size sensor and TV3050 is used with a 1/3" size sensor. **TV Lens** and its corresponding optical parameters are shown in Table 3.

Table 3 The order number and optical parameters of TV Lens

| Order number | Magnification | The maximum compatible sensor size |
|--------------|---------------|------------------------------------|
| TV3050 | 0.50X | 1/3" |
| TV3075 | 0.75X | 1/1.8" |
| TV3100 | 1.00X | 2/3" |

Note: Smaller image sensors can also be used for **TV Lens** with higher magnification, but in this way, the FOV of **MZO** will be reduced.

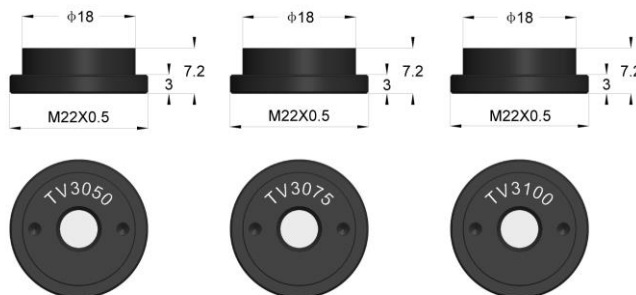


Figure 4 TV Lens with different magnification for TZM0880

1.3.4 Module Four: Bracket Adapter

Bracket Adapter (code **Add**) used for brackets with different mounting apertures and is sheathed in the middle of **MZO**. **dd** is the diameter of the bracket. Current available **Adds** include **A50** and **A45**. The installation method of the **Bracket Adapter** and bracket is shown in Figure 5.

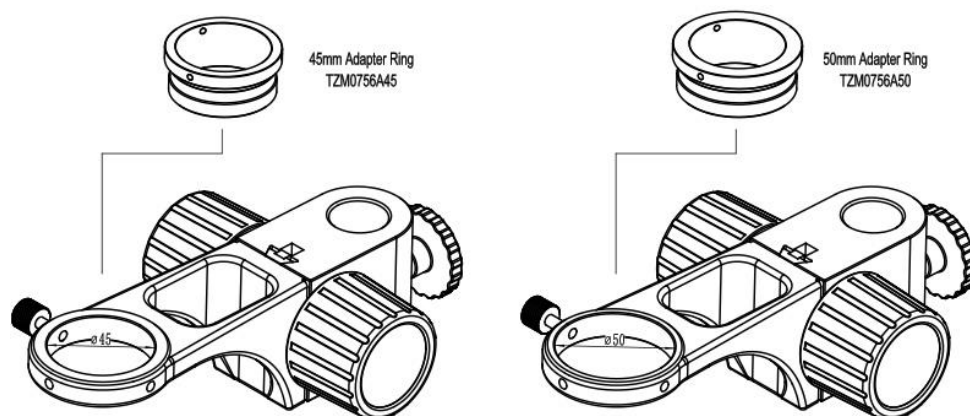


Figure 5 TZM0880 series' Bracket Adapter

1.3.5 Module Five: The Optional Module

The **Optional Modules** include **LED Direct Ring Light Module** (including **LED Direct Ring Light Module** and **LED Direct Ring Polarization Light Module**) and **Coaxial Light Module** (composed of **Coaxial Light Adapter** and **LED Spot Light**). The currently possible samples are shown in Table 4.

Table 4 The Optional Module: Light source module for TZM0880 series MZO

| Module | Order Number | Description |
|------------------------------|------------------|------------------------------------|
| LED Direct Ring Light Module | TZM0756DRL-65/85 | LED Direct Ring Light |
| | TZM0756DRPL-65 | LED Direct Ring Polarization Light |

| | | |
|---------------------------|---------------------|--|
| Coaxial Light Module | TZM0756CL+TZM0756SL | Coaxial Light Adapter + LED Spot Light(3W) |
| Transmissive Light Module | TZM0756TL | LED Transmissive Light |
| Power of Light Source | 40600014 | US:POWER-U-12V1A(Power Adapter American Standard) |
| | 40600015 | DE: POWER-E-12V1A(Power Adapter European Standard) |

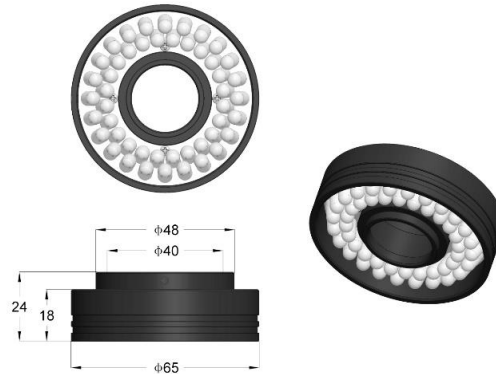


Figure 6 TZM0756DRL-65, the maximum diameter of LED reaches 65mm, and the diameter of 40mm hole in the center can be directly sleeved to the front end of the corresponding MZO to achieve uniform reflection lighting

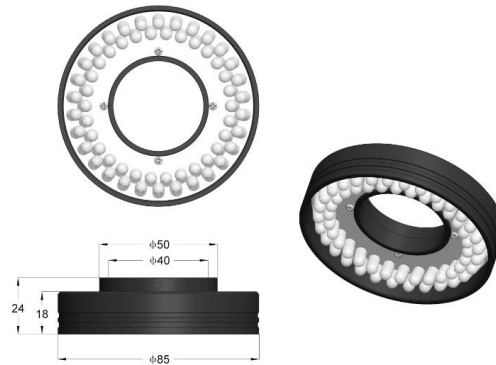


Figure 7 TZM0756DRL-85, the maximum diameter of LED reaches 85mm, and the diameter of 40mm hole in the center can be directly sleeved to the front end of the corresponding MZO to achieve uniform reflection lighting

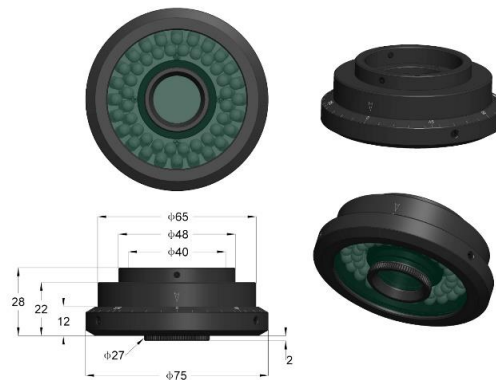


Figure 8 TZM0756DRPL-65, the maximum diameter of the LED reaches 65mm, and the small hole with a center diameter of 40mm can be directly sleeved on the front end of the corresponding MZO. By rotating the polarizing plate, uniform illumination of linear polarized light in different directions can be realized

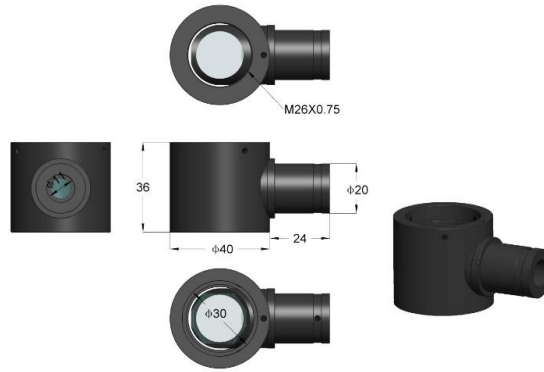


Figure 9 TZM0756CL Coaxial Light Adapter with 11mm hole for LED Spot Light

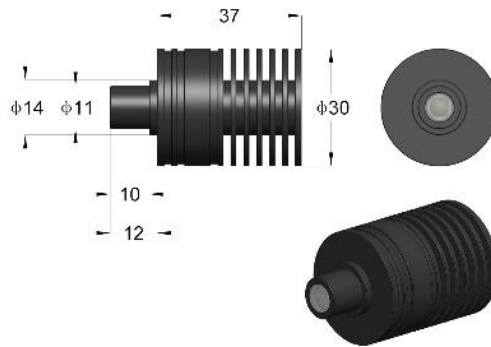


Figure 10 TZM0756SL, LED Spot Light can be directly connected with Coaxial Light adapter(TZM0756CL) to realize coaxial light illumination

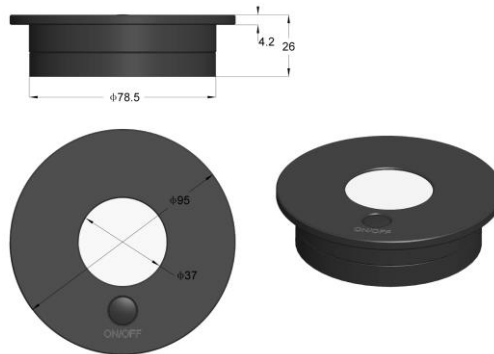


Figure 11 TZM0756TL LED Transmission Light source with 37mm illumination range

1.3.6 The Nomenclature of TZM0880 Series MZO

We have a special naming format for TZM0880 series MZO, we name TZM0880 series products with a series module code, which is a combination of code of different modules and connected with crossbar “-” to form a comprehensive naming. With this paradigm, user can understand the meaning and function briefly. For example, if a MZO is composed of Middle Zoom Module TZM0880D, Auxiliary Lens W3100(1.00x), TV Lens TV3050(0.5x) and 50mm Bracket Adapter A50, the final MZO can be expressed as:

TZM0880D-W3100-TV3050-A50

Currently available TZM0880a: TZM0880, TZM0880D (D represents Detent);

Currently available **W3bbb**: W3100, W3035, W3050, W3075, W3150;

Currently available **TV3ccc**: TV3100, TV3050, TV3075;

Currently available **Add**: A50 and A45;

Currently available Optional Modules are **LED Direct Ring Light Module** (including **LED Direct Ring Light Module** and **LED Direct Ring Polarization Light Module**) and **Coaxial Light Module** (composed of **Coaxial Light Adapter** and **LED Spot Light**).

The other **Auxiliary Lens** with **b.bb** magnification, **TV Lens** with **c.cc** magnification and **Bracket Adapter** with **dd mm** size can be customized according to the requirements.

1.3.7 The Configuration of TZM0880 with Five Modules

TZM0880 series **MZO** can be combined with **Camera Module** (optional) to form a digital monocular zoom microscope.

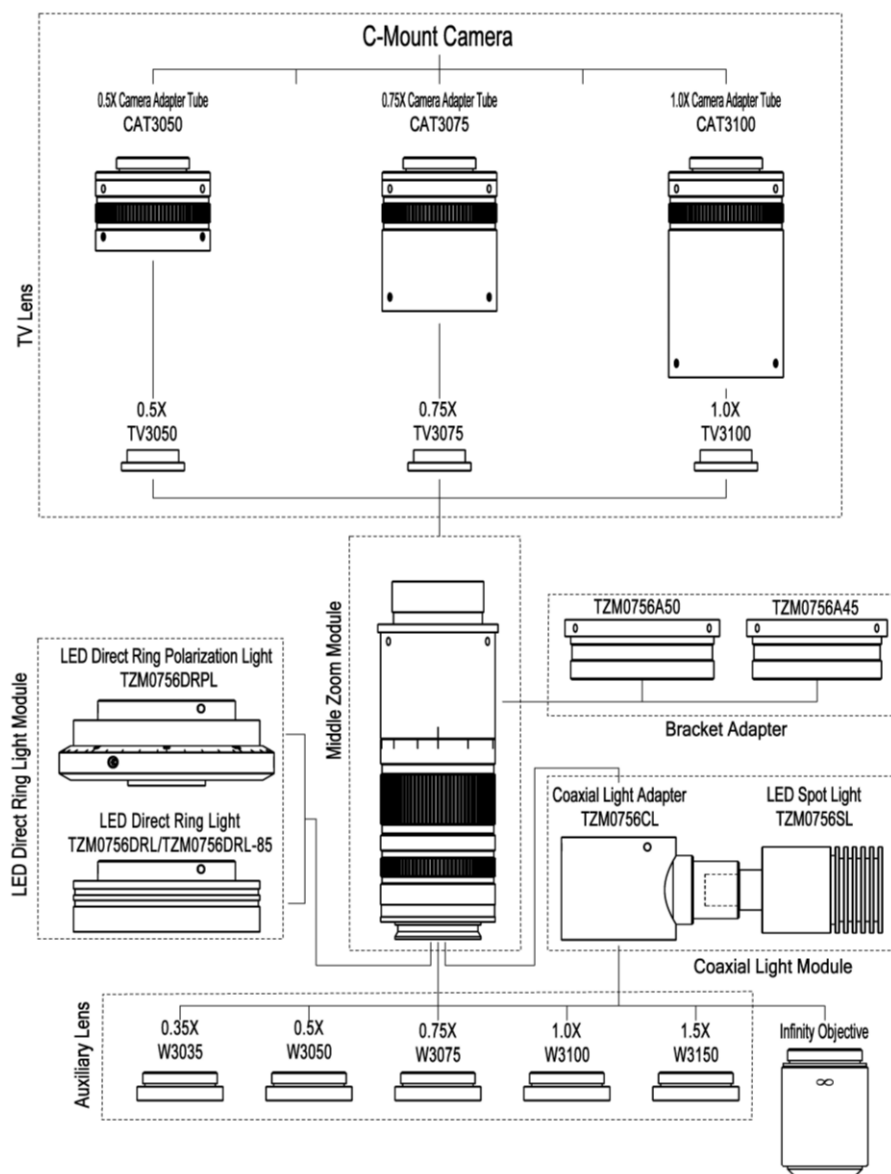


Figure 12 The Five Modules and TZM0880 MZO

1.4 The Basic Characteristics of TZM0880 Series MZO

The basic module of **TZM0880** series monocular zoom object is **TZM0880-W3100-TV3100-A50**. Its main features are shown below:

- With 0.8X~8.0X zoom range, 10X zoom ratio
- Larger NA: 0.02-0.11 (When using 1.00x **Auxiliary Lens**)
- Higher resolution: 15.3um-2.45um (When using 1.00x **Auxiliary Lens**)
- Larger field of view: 0.89mm-42.86mm (Object plane)
- Larger sensor size: 2/3" (When using 1.00x **TV Lens**)
- Working distance range: 56mm-248mm
- With adjustable BFL, parfocal in zoom range
- With adjustable center, the image center remains unchanged from 8.0x to 0.8x
- Compatible with infinity objectives (both biological and metallographic)
- Compact size: 194 mm (length) × 40 mm (diameter)
- **Auxiliary Lens** with 0.35x, 0.50x, 0.75x, 1.00x, 1.50x magnification (Optional)
- **TV Lens** with 0.50x, 0.75x, 1.00x magnification (Optional)
- Brightness adjustable **LED Direct Ring Light** (Optional)
- Brightness adjustable **LED Direct Ring Polarization Light** (Optional)
- Brightness adjustable **LED coaxial illumination** (Optional)
- **LED Transmissive Light** (Optional)
- 45mm or 50mm **Bracket Adapter** (Optional)

1.5 The Optical Specifications of TZM0880 with Different Auxiliary Lens and TV Lens

The optical specifications of **TZM0880** with different **Auxiliary Lens** and **TV Lens** are shown in Table 5. **Auxiliary Lens** and **TV Lens** with 1.00x are listed in the left-up cell. Its data is the basis of the other parameters in the whole table.

Table 5 TZM0880-W3100-TV3100 and its extensions

| Auxiliary Lens | Specification | TV3100 2/3"(D=11mm) | | TV3050 1/3"(D=6mm) | | TV3075 1/2"(D=8mm) | |
|---------------------------------|---------------|------------------------|-------|-----------------------|-------|-----------------------|-------|
| W3100 WD 86mm | PMAG | 0.80 | 8.00 | 0.40 | 4.00 | 0.60 | 6.00 |
| | DFOV/mm | 13.75 | 1.38 | 15.00 | 1.50 | 13.33 | 1.33 |
| | NA | 0.018 | 0.112 | 0.018 | 0.112 | 0.018 | 0.112 |
| W3035 WD 248mm | PMAG | 0.28 | 2.80 | 0.14 | 1.40 | 0.21 | 2.10 |
| | DFOV/mm | 39.29 | 3.93 | 42.86 | 4.29 | 38.10 | 3.81 |
| | NA | 0.006 | 0.039 | 0.006 | 0.039 | 0.006 | 0.039 |
| W3050 WD 174mm | PMAG | 0.40 | 4.00 | 0.20 | 2.00 | 0.30 | 3.00 |
| | DFOV/mm | 27.50 | 2.75 | 30.00 | 3.00 | 26.67 | 2.67 |
| | NA | 0.009 | 0.056 | 0.009 | 0.056 | 0.009 | 0.056 |
| W3075 WD 115mm | PMAG | 0.60 | 6.00 | 0.30 | 3.00 | 0.45 | 4.50 |
| | DFOV/mm | 18.33 | 1.83 | 20.00 | 2.00 | 17.78 | 1.78 |
| | NA | 0.014 | 0.084 | 0.014 | 0.084 | 0.014 | 0.084 |
| W3150 WD 56mm | PMAG | 1.20 | 12.00 | 0.60 | 6.00 | 0.90 | 9.00 |
| | DFOV/mm | 9.17 | 0.92 | 10.00 | 1.00 | 8.89 | 0.89 |
| | NA | 0.027 | 0.168 | 0.027 | 0.168 | 0.027 | 0.168 |

WD: Working Distance;

PMAG: Primary Magnification;

FOV: Field of View in the object side;

NA: Numerical Aperture;

Note: Infinity corrected objectives will limit system's usable zoom range due to uneven illumination.

1.6 The Dimension of TZM0880 Series MZO

The dimension of TZM0880 series MZO with different Auxiliary Lens and TV Lens are shown in Figure 13. The length of the MZO with different TV Lens will be slightly different. The length of TZM0880-W3100-TV3100-A50 is 194mm.

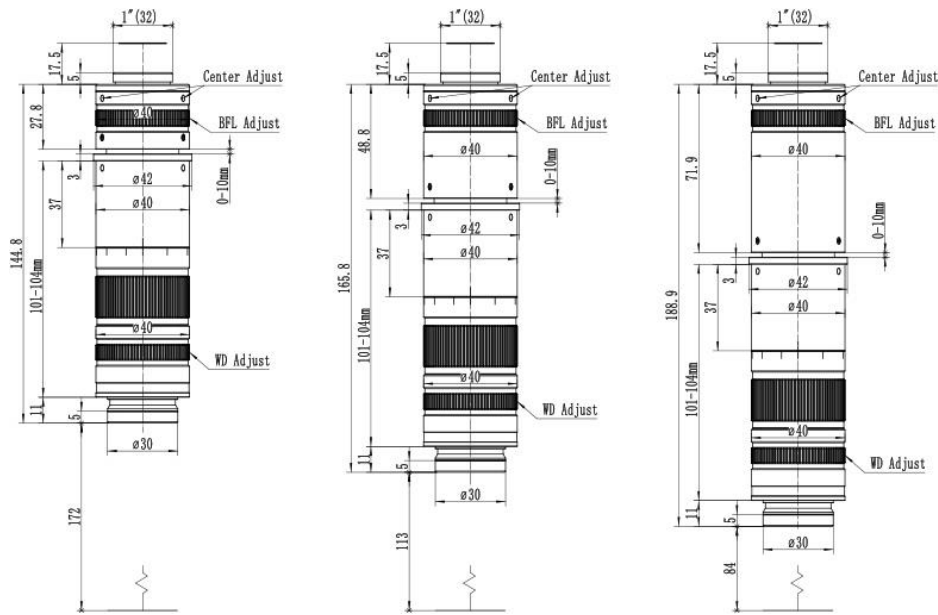


Figure 13 Dimension of TZM0880 series MZO with different Auxiliary Lens and TV Lens

a) TZM0880-W3050-TV3050; b) TZM0880-W3075-TV3075 c) TZM0880-W3100-TV3100

The dimensions of TZM0880-W3100-TV3100-A50 with different light are shown in Figure 14.

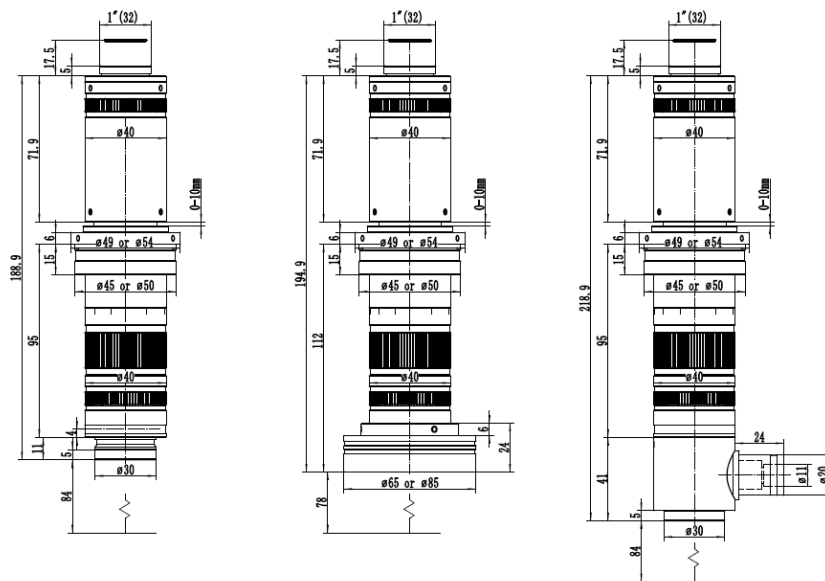


Figure 14 Dimensions of TZM0880-W3100-TV3100-A50 with different light module

a) TZM0880-W3100-TV3100-A50 without light module,

- b) TZM0880-W3100-TV3100-A50 equipped with Direct Ring Light Module,
c) TZM0880-W3100-TV3100-A50 equipped with Coaxial Light Module

1.7 How to Configure TZM0880 Series MZO

The corresponding parameters of TZM0880 series MZO are listed in Table 6. A specific combination can be determined according to the following steps.

1. Confirm the possible range of 1) FOV and 2) Working Distance in the object space to choose the Auxiliary Lens.
2. Choose the M26x0.705 to M20x0.705 Objective Adapter, if the M20x0.705 infinity objective is used.
3. Confirm the camera Image Area Size, it can be 1) Sensor Size (1/x in inch), 2) Image Diagonal Length, 3) Image Width or 4) Image Height to choose the right TV Lens.
4. Choose the 45mm or 50 mm adapter according to the diameter of the bracket hole diameter.
5. Choose the LED Direct Ring Light Module for the reflective illumination.
6. Choose the Coaxial Light Module if coaxial illumination is required.
7. Choose the Transmissive Light Module if transmissive illumination is required.
8. Choose the Camera Module.

We offer a variety of microscope cameras. Customers could get the detailed information of our cameras on the official website (<http://www.touptek.com/>) and choose the appropriate ones.

Table 6 The combinations of different the Auxiliary Lens, The Middle Zoom Module and the TV Lens. With this figure, user can configure Monocular Zoom Objective with different magnification to fit different sensor and different application

| Camera | Auxiliary Lens-Main Zoom Module-TV Lens | WD(mm) | PMAG | DFOV(mm) | NAO |
|---------|---|--------|------------|------------|-------------|
| 1/3inch | TZM0880-W3035-TV3050 | 248 | 0.14X-1.4X | 42.86-4.29 | 0.006-0.039 |
| | TZM0880-W3050-TV3050 | 174 | 0.2X-2.0X | 30.00-3.00 | 0.009-0.056 |
| | TZM0880-W3075-TV3050 | 115 | 0.3X-3.0X | 20.00-2.00 | 0.014-0.084 |
| | TZM0880-W3100-TV3050 | 86 | 0.4X-4.0X | 15.00-1.50 | 0.018-0.112 |
| | TZM0880-W3150-TV3050 | 56 | 0.6X-6.0X | 10.00-1.00 | 0.027-0.168 |
| 1/2inch | TZM0880-W3035-TV3075 | 248 | 0.21X-2.1X | 38.10-3.81 | 0.006-0.039 |
| | TZM0880-W3050-TV3075 | 174 | 0.3X-3.0X | 26.67-2.67 | 0.009-0.056 |
| | TZM0880-W3075-TV3075 | 115 | 0.45X-4.5X | 17.78-1.78 | 0.014-0.084 |
| | TZM0880-W3100-TV3075 | 86 | 0.6X-6.0X | 13.33-1.33 | 0.018-0.112 |
| | TZM0880-W3150-TV3075 | 56 | 0.9X-9.0X | 8.89-0.89 | 0.027-0.168 |
| 2/3inch | TZM0880-W3035-TV3100 | 248 | 0.28X-2.8X | 39.29-3.93 | 0.006-0.039 |
| | TZM0880-W3050-TV3100 | 174 | 0.4X-4.0X | 27.5-2.75 | 0.009-0.056 |
| | TZM0880-W3075-TV3100 | 115 | 0.6X-6.0X | 18.33-1.83 | 0.014-0.084 |
| | TZM0880-W3100-TV3100 | 86 | 0.8X-8.0X | 13.75-1.38 | 0.018-0.112 |
| | TZM0880-W3150-TV3100 | 56 | 1.2X-12X | 9.17-0.92 | 0.027-0.168 |

2 The Connection of TZM0880 Series MZO and Camera



Figure 15 TZM0880-W3100- TV3050+TZM0756DRL-65 (LED Direct Ring Light Module)+
HDMI Camera XCAMLite1080P A(Circular)+TPS-20A+SCRN

3 Pictures Taken with TZM0880 Series MZO

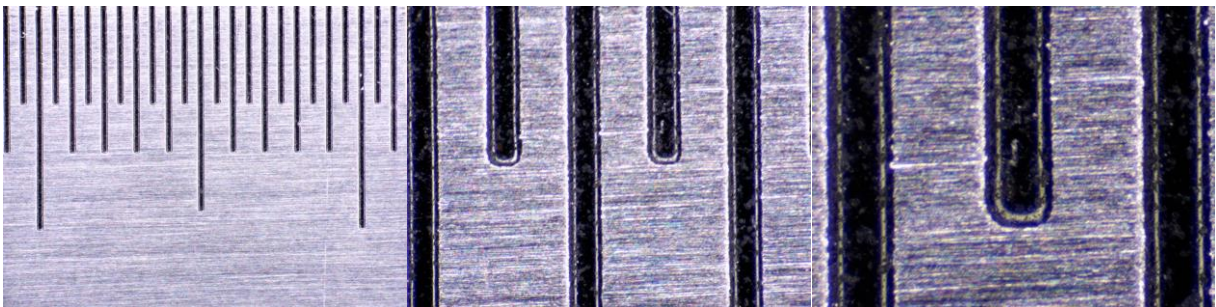


Figure 16 Pictures of steel ruler taken with TZM0880-W3100-TV3050 under LED Direct Ring Light illumination
left) 0.8X, middle) 4.0X, right)8.0X

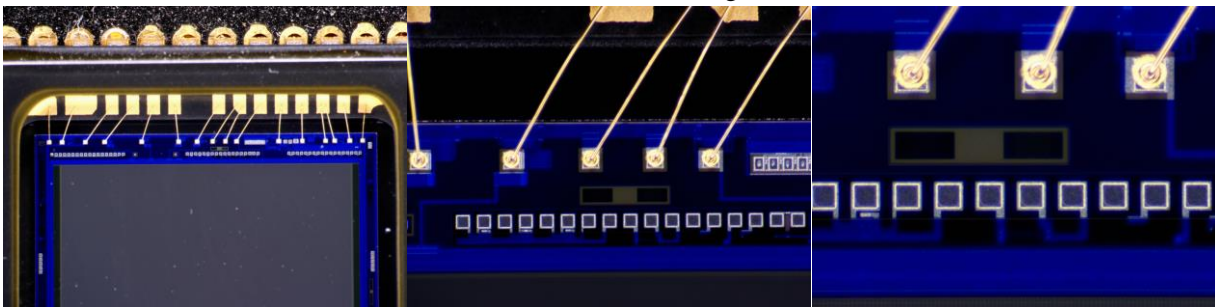


Figure 17 Pictures of CCD taken with TZM0880-W3100-TV3050
left) 0.8X, middle) 4.0X, right) 8.0X

4 TZM0880 Series MZO Packing List

The packing information of the [TZM0880](#) series MZO is as follows:



Figure 18 The TZM0880 Main Body, including Auxiliary Lens Module, Middle Zoom Module, TV Lens, Camera Adapter Tube and Bracket Adapter

The packing information of [TZM0756DRL-65/85](#) is as follows:



Figure 19 TZM0756DRL-65/85, including LED Direct Ring Light and Power Adapter

The packing information of [TZM0756DRPL-65](#) is as follows:



Figure 20 TZM0756DRPL-65, including LED Direct Ring Polarization Light and Power Adapter

The packing information of [Coaxial Light Module](#) is as follows:



Figure 21 TZM0880 Coaxial Light Module, including TZM0756CL (Coaxial Light Adapter), TZM0756SL (LED Spot Light) and Power Adapter

5 Order List of TZM0880 Series MZO

TZM0880 Series Monocular Zoom Object Package Order List

| Order number | | Quantity | Remarks |
|---|---------------------------|----------|---------|
| Package Name | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| | TZM0880□-W3□□□-TV3□□□-A□□ | | |
| Remarks | | | |
| Remarks: Users or salesmen can directly write the corresponding module name, number of sets, and other special requirements in the remarks with their own professional knowledge. | | | |

TZM0880 Order List of Other Accessories

| Module | Order number | Quantity | Remarks | Description |
|------------------------------|---------------------|----------|---------|--|
| Auxiliary Lens Module | W3035 | | | 0.35X Object Lens |
| | W3050 | | | 0.50X Object Lens |
| | W3075 | | | 0.75X Object Lens |
| | W3100 | | | 1.0X Object Lens |
| | W3150 | | | 1.5X Object Lens |
| | ON-XX | | | Biological Objective |
| | ON-YY | | | Metallographic Objective |
| | Objective Adapter | | | M26x0.706 to M20x0.706 |
| TV Lens | TV3050 | | | For 1/3" Sensor |
| | TV3075 | | | For 1/1.8" Sensor |
| | TV3100 | | | For 2/3" Sensor |
| Bracket Adapter | A45 | | | 45mm Bracket Adapter |
| | A50 | | | 50mm Bracket Adapter |
| | | | | |
| Coaxial Light Module | TZM0756CL+TZM0756SL | | | Coaxial Light Adapter + LED Spot Light(3W) |
| LED Direct Ring Light Module | TZM0756DRL-65 | | | LED Direct Ring Light (65mm outer diameter) |
| | TZM0756DRL-85 | | | LED Direct Ring Light (85mm outer diameter) |
| | TZM0756DRPL-65 | | | LED Direct Ring Polarization Light |
| Transmissive Light Module | TZM0756TL | | | LED Transmissive Light |
| Power of light source | 40600014 | | | POWER-U-12V1A(MSA-C1000IC12.0-12H-US), Power Adapter American Standard |
| | 40600015 | | | POWER-E-12V1A(MSA-C1000IC12.0-12W-DE), Power Adapter European standard |

TZM0880 Function Module Optional Order List

| Module | Order number | Quantity | Remarks | Description |
|---------------------------------|---------------------|----------|---------|---|
| Middle Zoom | TZM0880 | | | Ordinary Middle Zoom Module |
| Module | TZM0880D | | | Middle Zoom Module with detent |
| | | | | |
| Auxiliary Lens Module | W3035 | | | 0.35X Object Lens |
| | W3050 | | | 0.50X Object Lens |
| | W3075 | | | 0.75X Object Lens |
| | W3100 | | | 1.0X Object Lens |
| | W3150 | | | 1.5X Object Lens |
| | ON-XX | | | Biological Objective |
| | ON-YY | | | Metallographic Objective |
| | Objective Adapter | | | M26x0.706 to M20x0.706 |
| | | | | |
| TV Lens | TV3050 | | | For 1/3" Sensor |
| | TV3075 | | | For 1/1.8" Sensor |
| | TV3100 | | | For 2/3" Sensor |
| | | | | |
| Bracket Adapter | A45 | | | 45mm Bracket Adapter |
| | A50 | | | 50mm Bracket Adapter |
| Coaxial Light Module | TZM0756CL+TZM0756SL | | | Coaxial Light Adapter + LED Spot Light(3W) |
| LED Direct Ring Light Module | TZM0756DRL-65 | | | LED Direct Ring Light (65mm outer diameter) |
| | TZM0756DRL-85 | | | LED Direct Ring Light (85mm outer diameter) |
| | TZM0756DRPL-65 | | | LED Direct Ring Polarization Light |
| Transmissive Light Module | TZM0756TL | | | LED Transmissive Light |
| Power of Light Source | 40600014 | | | POWER-U-12V1A(American Standard) |
| | 40600015 | | | POWER-E-12V1A(European standard) |

6 Service

For customers who purchase MZO from local dealer, please contact your dealer for more information.

For technical support, please contact e-mail address: tphz@touptek.com.